

Analysis of Analgesics Using Atlantis T3

Waters Corporation



This is an Application Brief and does not contain a detailed Experimental section.

Abstract

This application brief highlights the analysis of analgesics.

Introduction

Compounds analysed in this application brief :

1. Acetaminophen
2. Caffeine
3. 2-Acetamidophenol
4. Acetanilide
5. Acetylsalicylic acid
6. Phenacetin

Experimental

Test Conditions

Column:	Atlantis T3, 4.6 x 150 mm, 5 µm
Part Number:	186003747
Mobile Phase A:	H ₂ O
Mobile Phase B:	ACN
Mobile Phase C:	1% HCOOH in H ₂ O, pH 2.3
Flow Rate:	1.0 mL/min
Injection Volume:	10 µL
Sample:	Acetaminophen (20 µg/mL), Caffeine (20

µg/mL), 2-Acetamidophenol (50 µg/mL),
Acetanilide (20 µg/mL), Acetylsalicylic acid
(50 µg/mL), Phenacetin (20 µg/mL) in H₂
O

Column Temperature: 30 °C

Sample Temperature: 15 °C

Detection: UV @ 260 nm

Sampling Rate: 5 point/sec


Filter Response: 0.2

Instrument: Waters Alliance 2695 with 2996 PDA

Gradient

Time (min)	Profile
	%A
0.00	75
10.00	30
11.00	75
15.00	75

Results and Discussion

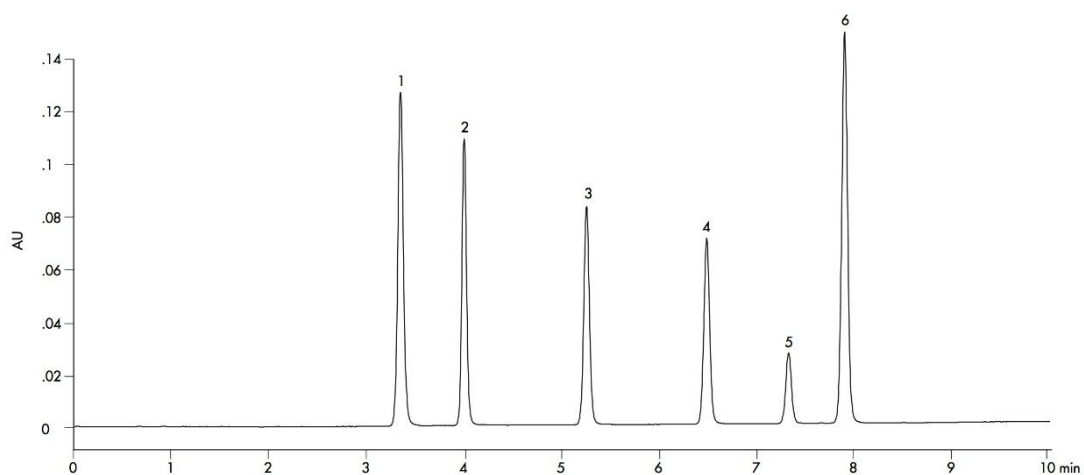

1. Acetaminophen, 2. Caffeine, 3. 2-Acetamidophenol, 4. Acetanilide, 5. Acetylsalicylic acid, 6. Phenacetin

Featured Products

Alliance HPLC System <<https://www.waters.com/534293>>

2998 Photodiode Array (PDA) Detector <<https://www.waters.com/1001362>>

WA60217, January 2009