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The screenshot displays the Peptide MAM software interface. On the left, a navigation menu includes 'Peptide MAM', 'Define Attributes', 'Define SSTs', and 'Set Processing Parameters'. The main area is divided into two sections: 'Peptide Identification' and 'New Peak Detection'.
Peptide Identification: This section allows users to set 'Mass Tolerance (ppm)' to 10 and 'Retention Time Tolerance (min)' to 0.5. Under 'Peak Processing', there are checkboxes for 'Ignore ions before' (set to 2.00 minutes) and 'Ignore ions after' (set to 52.00 minutes).
New Peak Detection: This section sets criteria for new peaks, including 'Fold change >=' (5.0), '%Base Peak Intensity >=' (0.1), 'Peptidic Isotope Profile Score >=' (75), and 'Retention Time Match Tolerance (min)' (1). There is also a checkbox for 'Exclude peptides with only +1 ions'.

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Reports

Report Content

Select the sections to include in your report:

- Injection List
- Blanks
- System Suitability
- Monitored Attributes
- New Peak Detection
- Acquisition Record

Generate Report

Peptide MAM Report

Generated By: waters_connect Administrator
 Generated Date: 29 Sep 2020 12:46
 Processing Date: 25 Sep 2020 13:06

Analysis
 Name: Peptide MAM
 Version: 2
 Location: Company\MAM\Demo data
 Description: bundle 34

Processing Method
 Name: Peptide MAM
 Version: 2
 Location: Company\MAM\Demo data
 Description: No Description

Injection List

#	Injection Name	Type	Acquisition Status
1	Blank_1	Blank	Complete
2	System suitability peptides_1_537	System Suitability	Complete
3	System suitability peptides_3_537	System Suitability	Complete
4	System suitability peptides_4_537	System Suitability	Complete
5	Blank_2	Blank	Complete
6	Control	Unknown	Complete
7	Unknown sample 1	Unknown	Complete
8	Unknown sample 2	Unknown	Complete
9	Unknown sample 3	Unknown	Complete
10	Blank_3	Blank	Complete
11	System suitability peptides_4_537	System Suitability	Complete
12	System suitability peptides_5_537	System Suitability	Complete
13	System suitability peptides_6_537	System Suitability	Complete
14	Blank_4	Blank	Complete

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Results - % Modified

Injection Name	DIQMPSPPTPLAS NPD Response S/N Ratio	DIQMPSPPTPLAS NPD Variation S/N Ratio	DIQMPSPPTPLAS NPD Response S/N Ratio	DIQMPSPPTPLAS NPD Variation S/N Ratio	DIQMPSPPTPLAS NPD Response S/N Ratio	DIQMPSPPTPLAS NPD Variation S/N Ratio
Control	1.26	0.57	0.62	1.47	0.22	0.22
Unknown sample 1	1.18	1.18	0.68	4.18	42.33	42.33
Unknown sample 2	1.24	0.85	0.82	1.36	42.76	42.76
Unknown sample 3	1.25	1.49	0.68	4.54	42.44	42.44

New Peak Detection

New Peak Detection Criteria

Criteria	Value
Peak Change	>1%
Baseline	>100ms
Isolated Intensity	>60000
Marking Tolerance (sigma)	1
Exclude peaks with only 1 event	Yes

Processing Parameters

Parameter	Value
System suitability criteria	2.00
Injection delay (min)	0.00

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ACQUITY HSS Detector

Serial Number: 15A042

Instrument Setup

Method: Report
 Location: Method 2005
 Last performed: Mar 10, 2020 10:16:41 EDT
 Location type: System

LC-MS Binary Solvent Manager

Software version: 3.1.2
 Checksum: 40268414
 Mean Argon gas pressure: 0.48 psi
 Mean Argon gas pressure: 0.47 psi
 Mean Argon gas pressure: 0.48 psi
 Mean Argon gas pressure: 0.48 psi

Sample Manager 71N

Software version: 3.1.2
 Checksum: 40268414
 Sample Name: 15A042
 Column Name: 15A042
 Column Station: 10
 Min Sample Temperature: 6.0°C
 Min Column Temperature: 6.0°C
 Min Column Temperature: 6.0°C
 Min Column Temperature: 6.0°C
 Auto Inhibit Enabled: Yes

UV Detector

Software version: 3.1.2
 Checksum: 40268414
 Wavelength: 214 nm
 Flow Cell Part Number: 010
 Flow Cell Part Number: 010
 Flow Cell Part Number: 010
 Flow Cell Part Number: 010
 Lamp On: 15A042
 Message:

System suitability peptides_2

Experimental Record

UV Detector

Wavelength mode: Single wavelength
 Wavelength: 214 nm
 Sampling Rate: 20 points/min

Channel A

Channel Wavelength: 214 nm
 Flow Flow Control Mode: Report
 Auto Zero On Wavelength Change: Maintain baseline
 Channel:

Acquisition

Using Report 1
 Data Mode: Acquisition
 Priority: Acquisition 1
 Disable Chem Warn: Yes

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1. Xu, Weichen *et al.* A Quadrupole Dalton-based Multi-Attribute Method for Product Characterization, Process Development, and Quality Control of Therapeutic Proteins. *mAbs* vol. 9,7 (2017): 1186–1196. doi:10.1080/19420862.2017.1364326.
2. Alt N, Zhang TY, Motchnik P, Taticek R, Quarmby V, Schlothauer T, Beck H, Emrich T, Harris RJ. Determination of Critical Quality Attributes for Monoclonal Antibodies using Quality by Design Principles. *Biologicals*. 2016 ;44:291–305. doi:10.1016/j.biologicals.2016.06.005. PMID:27461239.
3. Finkler C, Krummen L. Introduction to the Application of QbD Principles for the Development of Monoclonal Antibodies. *Biologicals*. 2016;44:282–90. doi:10.1016/j.biologicals.2016.07.004. PMID:27546184.
4. Rogers, Richard S *et al.* Development of a Quantitative Mass Spectrometry Multi-Attribute Method for Characterization, Quality Control Testing and Disposition of Biologics. *mAbs* vol. 7,5 (2015): 881–90. doi:10.1080/19420862.2015.1069454.
5. Li, W, Kerwin, J. L, Schiel, J, Formolo, T, Davis, D, Mahan, A, Benchaar, S. A. State-of-the-Art and Emerging Technologies for Therapeutic Monoclonal Antibody Characterization Volume 2. *Biopharmaceutical Characterization: The NISTmAb Case Study*. January 1, 2015 , 119–183.

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© 2023 Waters Corporation. All Rights Reserved.

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